RR #2, Warton, Ontario, CANADA N0H 2T0

Toll Free (Canada / USA)
Telephone
Facsimile

800-567-3556
519-534-1080
519-534-1088

□ BDC2002

□ BDC201

ENGLISH

Instruction Manual
pages 2-17

ESPAÑOL

Manual de instrucciones
páginas 34-49

FRANÇAIS

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DEUTSCH

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Garantie

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CONTENTS

| | |
|---|-------|
| Safety Instructions..... | 3 |
| General Instructions..... | 4 |
| Packing List..... | 5 |
| BDC2002 Description..... | 6 |
| BDC2010 Description..... | 7 |
| Technical Data BDC2002..... | 8 |
| Technical Data BDC2010..... | 9 |
| Installation in the Laboratory..... | 10 |
| Environmental Operating Conditions..... | 10 |
| BDC2002 General Operating Instructions..... | 11-12 |
| BDC2010 General Operating Instructions..... | 13-15 |
| Routine Cleaning, Service, Transport & Storage..... | 16 |
| Warranty and Product Liability..... | 17 |

SAFETY INSTRUCTIONS

Each operator should read all instruction before using this device.

The stirrer must be securely fixed to a stable support. Mount to a Caframo stand using a Caframo heavy duty clamp. If other stands or clamps are used, the unit must be secured such that it will not fall if tipped.

CAUTION : This is not an explosion proof stirrer. Do not use in a highly flammable or explosive environment.

Spinning paddles or impellers can cause severe personal injuries. Operators must use extreme care and good judgment when mixing at any speed.

All mixing paddles and impellers must be in good condition with straight shafts. If the stirrer vibrates at high speeds check the paddle shaft for damage and repair or replace if needed.

Extreme care must be taken when mixing chemicals to ensure that no chemicals are splashed outside the mixing vessel. Care must be taken when changing to faster mixing speeds.

Ensure that the mixing impeller does not contact the containment vessel.

Do not operate while standing in water. Keep the unit dry and do not immerse any part, except the mixing paddle into any liquids. Protect from splashing.

Ensure that no loose clothing, jewelry, or hair can become entangled in any rotating parts. A fast spinning chuck can cause injury to operator.

Power can be interrupted to the stirrer by pressing the power button. In an emergency, disconnect the mains cord.

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GENERAL INSTRUCTIONS

Read safety instructions.

Unpack the stirrer carefully. Ensure the stirrer is not damaged and verify the contents against the packing list. If the stirrer is damaged or the contents do not correspond with the packing list, notify your supplier immediately.

Read the manual carefully! Make sure that all users read this manual thoroughly before operating the stirrer. For further questions or inquiries, please contact:

Caframo Ltd.

R.R. #2, Warton, Ontario, Canada N0H 2T0
Tel : (519) 534-1080 Fax : (519) 534-1088

Toll free in USA and Canada :

Tel : (800) 567-3556 Fax : (800) 209-6786

Outside North America :

Email : labsales@caframo.com

Web Site :

www.caframo.com

APPROVALS

North American

CAN/CSA-C22.2 No. 1010-1:1993 (R1999)
ANSI/ISA S82.01-1994

European

EN61326 (EN55011, conducted and radiated emissions, EN61000-3-2, EN61000-4-2, EN61000-4-3, EN61000-4-4, EN61000-4-5, EN61000-4-6)

CSA
ANSI
ISA
CE
ITS
ETL

Canadian Standards Association
American National Standards Institute
Instrument Society of America
Communauté Européenne (European Community)
Intertek Testing Services (www.worldlab.com)
Electrical Testing Laboratory

PACKING LIST

Compare the contents of the shipment to the items on the following list to ensure all parts are received with the stirrer. Do not discard the container and packing materials until all parts are accounted for. It is recommended that the original packing materials be kept for transportation and storage.

- G CompactDigitalJ BDC2002 Stirrer with chuck or
- G ReversingDigitalJ BDC2010 Stirrer with chuck

- G Power Supply with 115 volt cord end or
- G Power Supply with 220 volt cord end

- G Support Bar (including allen screw, lock washer and key)

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TECHNICAL DATA BDC2002

Technical specifications subject to change without notice.

| Model | BDC2002 |
|--|---|
| Speed Range | 40 - 2002 rpm |
| LED Display Accuracy | " 1 rpm or 3% of reading, 4 digit |
| Maximum Torque | 70 N-cm, (100 in-oz) |
| Input to Power Supply | AC 100-240 volts, 50/60 Hz, 125-175 VA (Must use Caframo supplied power supply) |
| Electrical Output from Power Supply to Stirrer | DC 24 V, 2.7 A |
| Mechanical Output Power | 1/15 hp, 50 watts, (Brushless DC Motor) |
| Weight Stirrer Power Supply | 3.0 kg (7 lbs) 0.5 kg (1 lb) |
| Maximum Volume | 25 litres (6.6 US gallons, 5.3 Imp Gallons) |
| Maximum Viscosity | 15000 mPa.s (15000 cps) -Alike sour cream@ |
| Molded Front Housing | Flame retardant, chemical resistant, high impact nylon. Recyclable (PA). |
| Cast Back Housing | Epoxy coated zinc |
| Combined Housing | Splash Proof |
| Key Pad | Waterproof, chemical resistant polyester |
| StirlightJ | 5 Ultra Bright LED=s, 100,000 hour life |
| Power Cord Lengths | 193 cm (6') from Stirrer to Power Supply-DC 193 cm (6') from Power Supply to wall-AC |
| Adjustable Chuck | Fits onto 13mm (2 inch) Ø output shaft of stirrer. Adjusts up to 10 mm (3/8"). |

| Model | BDC2010 |
|--|---|
| Speed Range | 40 - 2010 rpm |
| Functions | Countdown Timer Clockwise Counter Clockwise xRx Agitation |
| LED Display Accuracy | " 1 rpm or 3% of reading, 4 digit |
| Maximum Torque | 100 N-cm, (142 in-oz) |
| Input to Power Supply | AC 100-240 volts, 50/60 Hz, 2.5 A (Must use Caframo supplied power supply) |
| Electrical Output from Power Supply to Stirrer | DC 24 V, 4.75 A, 90 W switching power supply |
| Mechanical Output Power | 1/10 hp, 70 watts, (Brushless DC Motor) |
| Weight | Stirrer 3.0 kg (7 lbs) Power Supply 0.6 kg (1.2 lbs) |
| Maximum Volume | 25 litres (6.6 US gallons, 5.3 Imp Gallons) |
| Maximum Viscosity | 20,000 mPa.s (20,000 cps) -Alike chocolate syrup@ |
| Molded Front Housing | Flame retardant, chemical resistant, high impact nylon. Recyclable (PA). |
| Cast Back Housing | Epoxy coated zinc |
| Combined Housing | Splash Proof |
| Key Pad | Waterproof, chemical resistant polyester |
| StirlightJ | 5 Ultra Bright LED=s, 100,000 hour life |
| Power Cord Lengths | 193 cm (6') from Stirrer to Power Supply-DC 193 cm (6') from Power Supply to wall-AC |
| Adjustable Chuck | Fits onto 13 mm (2 inch) Ø output shaft of stirrer. Adjusts up to 10 mm (3/8"). |

Technical specifications subject to change without notice.

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INSTALLATION IN THE LABORATORY

Installing the support bar

Remove the allen screw and lock washer from the support bar. Insert the support bar into the back of the stirrer so that the holes line up for the allen screw. Secure with allen screw and lock washer and tighten with allen key.

Clamp the stirrer to a Caframo stand on a stable, level surface. See safety instructions on page 3.

Connect stirrer to supplied Caframo power supply. Connect power supply to a grounded electrical outlet.

Impeller attachment

Insert the impeller shaft into the open chuck and adjust the impeller to the desired height. The hollow shaft allows the impeller shaft to come through the top of the stirrer if necessary.

CAUTION : Operating a freely rotating impeller shaft through the top of the stirrer is not safe. Do not allow the impeller shaft to touch or extend through the rubber gasket while the motor is powered on. The gasket is for sealing purposes only. Tighten the chuck's outer ring by hand and then secure with the chuck key.

ENVIRONMENTAL OPERATING CONDITIONS

The BDC2002 and BDC2010 stirrers must operate in the following conditions :

Indoors

1. Altitudes up to 6500 ft. (2000 m).
2. Temperatures from 41°F to 104°F (5°C to 40°C)
3. Maximum relative humidity 80% for temperatures up to 88°F (31°C) decreasing linearly to 50% relative humidity at 104°F (40°C).
4. Pollution degree 2 in accordance with IEC 664
"Normally only non conductive pollution occurs. Occasionally, however, temporary conductivity caused by condensation must be expected."

BDC2002 GENERAL OPERATING INSTRUCTIONS

Make sure the stirrer is securely mounted as specified in the installation instructions. Ensure the mixing paddle or impeller is securely attached to the desired position.

The stirrer is now ready for use to mix liquids and liquid/solid solutions. Appropriate impeller and container specifications must be determined by the operator, ensuring that all safety instructions are followed.

Button Operating Functions

Power Button - Turns the stirrer motor on and off.

Increase Button - Increases the speed by one rpm. If pressed within five seconds of the stirrer being turned on it will stir at the last speed before powered off.

Decrease Button - Decreases the speed by one rpm. If pressed within five seconds of the stirrer being turned on it will stir at 40 rpm.

Light Power Button - Turns the light on and off. The stirrer must be plugged in for the light to function.

Safety is built into this stirrer. When it is turned on it pauses for 5 seconds before it starts to mix. The last speed used is displayed on the LED. The user has three options.

1. Do nothing. The stirrer will start to stir at 40 rpm after five seconds.
2. Press the  and the stirrer will start to stir at the last speed used before powered off. This speed is displayed.
3. Press the  and the stirrer will start to stir at 40 rpm.

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LED Display Displays desired rotating speed in rpm. A “soft start” has been incorporated into the stirrer. When the increase or decrease buttons are pressed and held, the speed ramps smoothly so that getting to the speed you require is quick and easy.

Display Codes

- OL** Indicates an overload condition. The load is too great for the stirrer and must be reduced. To reset stirrer turn power off and back on.
- OH** Indicates an overheat condition. The operating conditions have caused the stirrer to overheat. Reduce stirrer loads and / or environmental temperature to continue mixing. To reset stirrer turn power off, allow stirrer to cool and then turn back on.

BDC2010 GENERAL OPERATING INSTRUCTIONS

Make sure the stirrer is securely mounted as specified in the installation instructions. Ensure the mixing paddle or impeller is securely attached to the desired position.

The stirrer is now ready for use to mix liquids and liquid/solid solutions. Appropriate impeller and container specifications must be determined by the operator, ensuring that all safety instructions are followed.

Keypad Functions

Power Button - Turns the stirrer on and off.

Multi-Function Button – Sets three parameters - speed (rpm), countdown timer (minutes) and xRx agitation (seconds)

Increase Button - Increases the speed by one rpm or the countdown timer by one minute or the xRx agitation by one second.

Decrease Button - Decreases the speed by one rpm or countdown timer by one minute or the xRx agitation by one second.

Reverse – Reverses the direction that the impeller is turning. The impeller turns clockwise by default.

Start/Pause Button – Allows the stirrer to be paused and started again. Use this button to start the stirrer when ready.

Light Power Button - Turns the light on and off. The stirrer must be plugged in for the light to function.

Four Digit LED Display – Displays either the speed (rpm), countdown timer (minutes) or xRx agitation (seconds). The four decimal points are used to indicate the following parameters. These do not act as numeric decimal points.

â indicator that speed is displayed

â indicator that speed is displayed

ã indicator that time is displayed

ã indicator that time is displayed

ä indicator that cycle time is displayed

ä indicator that cycle time is displayed

↻ indicator that stirrer is turning counter clockwise

- During set-up the indicator light is solid. When the stirrer is running, the indicator light flashes.

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- G Spannfutter

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Clockwise Stirring

Turn the stirrer on by pressing the power button. The display will read 40 rpm and the speed indicator light will be illuminated. Increase to the desired speed and press the start/pause button. The stirrer will start turning in a clockwise direction. The speed indicator light in the display will flash and the set speed in rpm will be displayed.

Counter Clockwise Stirring

Set up by following the clockwise stirring instructions above. Press the reverse button. The reverse button can be pressed while the stirrer is running to start the stirrer in a counter clockwise direction. The indicator light above the reversing icon will be illuminated when the stirrer is turning in a counter clockwise direction.

Countdown Timer

Turn the stirrer on by pressing the power button. Set the desired speed by pressing the “+” button. Press the multifunction button until the indicator light for time is illuminated. A “0” will appear in the display. Press the “+” to set the countdown timer in minutes. Press the start/pause button to start. Stirring will start. Time or speed can be displayed. If time is selected, the minutes remaining are displayed. Press the multifunction button to display speed. To stir in the opposite direction, press the reverse button. Stirring stops when time is completed. Three beeps will sound.

xRx Agitation Cycle

Turn the stirrer on by pressing the power button. Set speed as above and countdown timer if desired. Press the multifunction button once more until the indicator light for cycle is illuminated. A “0” will appear in the display. Press the “+” to set the number of seconds the impeller will rotate in one direction before changing directions. Press the start/pause button.

For example: If set for 5 seconds the stirrer impeller will turn clockwise for 5 seconds at the set speed then stop and change to turn counter clockwise for 5 seconds. This mixing pattern will continue until the pause button is pressed or until the time is completed (if set).

The display will show the cycle seconds counting down during each direction unless speed or the countdown timer is selected.

When the stirrer is powered on, the default speed is 40 rpm, countdown timer is 0 minutes, and the xRx agitation cycle is 0 seconds.

If the stirrer is kept powered on it will remember the original values set for speed, the countdown timer and the xRx agitation cycle so that when the start/pause button is pressed again the operations will repeat.

Speed can be changed (increased or decreased) while the stirrer is running. Press the multifunction button so the speed is displayed. Increase or decrease speed by pressing the "+" or "-" buttons.

The countdown timer and the xRx agitation cycle can be changed only when the stirrer is stopped and the indicator lights are solid. If the indicator lights are flashing the parameters cannot be changed. To reset the stirrer, power the stirrer off and then on again.

The multifunction button can be pressed when the stirrer is running to check the status of each function.

For safety, a "soft start" has been incorporated into the stirrer. The stirrer will ramp up to reach the set speed to minimize splashing.

In the event of a power outage, the stirrer will stop and set parameters will go back to their default values.

Display Codes

- OL Indicates an overload condition. The load is too great for the stirrer and must be reduced. Press the power button and enter in new parameters and start.

- OH Indicates an overheat condition. The operating conditions have caused the stirrer to overheat. Reduce stirrer loads and/or environmental temperature to continue mixing. To reset stirrer turn power off, allow stirrer to cool and then turn back on.

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ROUTINE CLEANING, SERVICE, TRANSPORT & STORAGE

Cleaning

Using a mild soap and water solution, wipe the housing and control panel of the stirrer with a damp cloth.

Note :

Do not use chlorine bleach, chlorine-based cleanser, abrasives, ammonia, steel wool or scouring pads with metal content or similar harsh solvents or abrasives. These may damage the surface of the instrument.

Service

Do not attempt to service or repair a Caframo stirrer. **If the stirrer housing is opened the warranty becomes void.** Contact Caframo for return authorization and return instructions. See page 4.

Transport and Storage

1. Clean stirrer so that it is free from any materials which may be harmful to the health. Provide a MSDS or other safety data sheets for all chemicals used in conjunction with this stirrer where appropriate.
2. Place the unit and its parts into the original packing or a container with the necessary protection to prevent damage during transport. Seal the original packing or container with packing tape.
3. Store the packed unit in a dry place.

CAUTION :

Failure to clean, service, and handle the stirrer as outlined can cause it to fail which could impair the safety of the user.

WARRANTY AND PRODUCT LIABILITY

Warranty

Caframo warrants the herein described product of its manufacture for three years from receipt of the equipment by the purchaser, against defects in material and workmanship. Warranty registration is available on-line. Point your browser to (www.caframo.com/warranty.htm) or fill out the supplied warranty registration card.

This limited warranty covers parts and labor.

If the stirrer housing is opened the warranty becomes void.

This warranty may only be altered by a specifically published amendment. No individual has authorization to alter the provisions of this warranty policy or its amendments. Expendable items are not covered by this warranty. Damages due to corrosion or accidental breakage are not covered.

Product Liability

Under no circumstances shall Caframo be liable for indirect, consequential or special damages of any kind.

Caframo reserves the right to make technical changes without prior notice.

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